

HPQA™, Helipath Quick Action Stand Firmware Update


Steps to update firmware

1. Insert a USB Flash Drive with the desired firmware version into the USB A Port on the back of the HPQA.




Be sure that no other HPQA firmware versions are on the flash drive.

2. From the Main Menu, select the Settings Icon

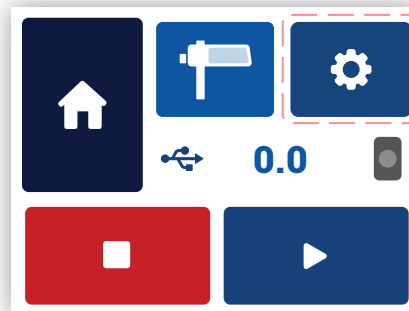

Settings Icon

3. Press the Firmware Update button.

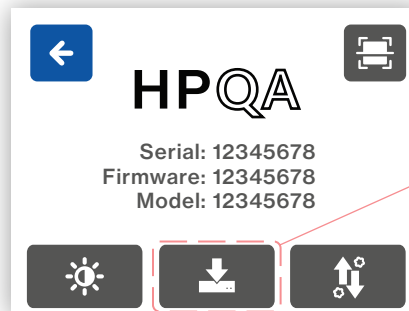

Firmware Update Button

4. The current version of firmware will be shown on the screen.
5. Update the firmware by pressing the circular download button on the firmware update screen.



Download Button

6. Wait for the HPQA to finish updating before pressing any buttons or moving the Joystick.



Firmware update may take a couple minutes. **DO NOT TURN THE HPQA OFF.**

7. From the Main Menu, select the Settings Icon and confirm Firmware version.